

### Agilent 5977A Series GC/MSD System

**Data Sheet** 



### 5977A GC/MSD Overview

The Agilent 5977A Series GC/MSD builds on a 45-year tradition of leadership and innovation, bringing together the technologies of the industry's best GC and MS systems. This advanced instrument provides a higher plane of productivity and confidence, including increased sensitivity, superior workflow, and software tools that simplify method optimization and lower operating costs.

# The Agilent GC/MSD System features Agilent 7890B GC, the most reliable and versatile gas chromatograph

- Precise pneumatic and temperature control for reproducible chromatographic performance.
- Low dead volume inert Capillary Flow Technology devices with various configurations for more productivity and more functionality.
- Sample Introduction systems to cover the widest range of samples for GC/MS analysis.

## Agilent 5977A GC/MSD, the most sensitive and robust mass selective detector

- · Extractor El ion source for high sensitivity, 10fg OFN IDL.
- Heated monolithic quartz gold quadrupole (heatable up to 200 °C) for rapid elimination of contamination to keep the analyzer clean.
- · Triple Axis Detector (TAD) for eliminating neutral noise.
- Scan Speed up to 20,000 u/sec (Extractor Ion Source).
- · Optional dry pump for oil-free operation.
- · Agilent TwisTorr 304 FS turbo pump optimized for helium and hydrogen.



#### Versatile workflow tools

- MassHunter Acquisition software with similar look and feel of MSD Productivity ChemStation Acquisition for efficient creation, update and use of GC/MS methods.
- Each system is delivered with both MassHunter Data Analysis and classic MSD ChemStation Data Analysis.

#### **GC-MS** direct communication

 Improved direct communication between the Agilent 7890B Series GC and 5977A MSD for optimized productivity and system protection.

### **Eco-friendly GC/MSD**

 The automated Sleep feature for reduction of electrical power and gas consumption with the Agilent DS42 inverter roughing pump, Agilent turbo pump and GC-MS direct communication.

### **Complete inert pathway**

 Prevents analyte breakdown from injection through detection for greater reproducibility and confidence in results.

### **Complete application solutions**

 Agilent analyzers and application kits for generating results immediately after installation.

#### **10 Year Value Promise**

 Guaranteed support for 10 years from the date of purchase, or Agilent will provide credit for the residual value of the system toward a model upgrade.

### **Agilent 5977A Series MSD System Specifications**

#### Mass spectrometer

| Mass spectrometer  |   |
|--|---|
| Ionization mode (standard)                                   | EI  |
| Ionization modes (optional)                                  | PCI, NCI, and EI acquisition with CI source   |
| El ion source type   | Noncoated Inert Extractor El source<br>Noncoated Inert El source  |
| Electron energy  | 5–241.5 eV  |
| Emission current   | 0–315 μΑ  |
| CI gases   | Dual gas inlet  |
| Filaments  | Dual for EI, single for CI  |
| Transfer line temperature                                    | 100–350 °C  |
| Ion source temperature                                       | 150–350 °C  |
| Quadrupole temperature                                       | 106–200 °C  |
| Mass filter  | Heated monolithic hyperbolic quadrupole   |
| Ion Source - Mass filter interface                           | Dynamically rampable entrance lens  |
| Mass range   | 1.6 to 1050 u   |
| Mass resolution  | Unit mass   |
| Mass axis stability  | Better than 0.10 u/48 hrs   |
| Detector   | Triple Axis Detector with long life triple channel EM   |
| Dynamic range (electronic)                                   | 10 <sup>6</sup>   |
| Scan rate <sup>1</sup>                                       | Up to 20,000 u/s (Extraction tune only), 12,500u/sec  |
| SIM  | 60 ions × 100 groups  |
| Pumping system (Helium)                                      | 65 L/s Diffusion Pump and 255 L/s Turbo Pump with 2.5 m <sup>3</sup> /hr mechanical pump  |
| Maximum recommended analytical gas flow                      | 1.5 mL/min (Diffusion Pump) 4 mL/min (Turbo Pump)   |
| Recommended maximum sustained column flow to MS <sup>2</sup> | 3 mL/min (Diffusion Pump) 50 mL/min (Turbo Pump)  |
| Instrument control   | MassHunter data system and local control panel  |
| Maintenance accessibility                                    | Source (including filaments, lenses), mass filter <sup>3</sup> , and detector on removable plate for easy access                                    |
| Maintenance scheduling                                       | Early Maintenance Feedback (EMF) for GC and MSD, user-defined maintenance schedule, with display of current status                                  |
| Power save<br>(Turbo Pump model only)                        | Sleep mode available for Turbo Pump and DS42i Executable Sleep/Wake from the panel of GC, Menu item in the MassHunter Software and Sequence Keyword |
| Support life   | Guaranteed support for 10 years   |
|  |   |

<sup>&</sup>lt;sup>1</sup>As scan rate increases, sensitivity will decrease and resolution may degrade.

<sup>&</sup>lt;sup>2</sup>High flow rate into a fixed ion source will cause a loss in sensitivity

<sup>&</sup>lt;sup>3</sup>The heated quadrupole mass filter should not require maintenance, but if maintenance is required, it should be performed by an Agilent service engineer

### Gas Chromatograph (Agilent 7890B GC)

| Autosampler                        | Agilent 7693, 7650, CombiPAL, 7697 and many third party autosamplers   |
|------------------------------------|--|
| Liner replacement                  | Compatible with turn-top system standard with 7890B GC   |
| Inlet (up to 2)                    | Split-splitless inlet (standard), Multimode Inlet (MMI), Cool On<br>Column (COC) inlet and others (optional) |
| Oven temperature                   | Ambient +4 – 450 °C  |
| Oven ramps/plateaus                | 20/21 with negative ramps allowed  |
| Carrier gases                      | Helium and hydrogen  |
| Electronic pneumatic control (EPC) | EPC split/splitless, septum purge and auxiliary EPC modules  |
| Carrier gas control modes          | Constant flow, constant pressure, programmable flow and programmable pressure                                |
| Pressure range                     | 0–100 psi (standard), 0–150 psi (optional) with 0.001 psi resolution   |
| Flow control                       | Compatible with optional capillary flow device controller  |

#### **Data System**

| •   |  |
|---|--|
| Acquisition Software  | MassHunter GC/MS Acquisition   |
| Simultaneous MS and GC detectors                                    | Up to two MS (SIM/scan) and four GC detector signals   |
| SIM/Scan  | Automated SIM setup and synchronous SIM/Scan operation   |
| Chemical ionization setup   | Electronic mass flow control of reagent gases  |
| High-mass confirmation  | Verification test kit (optional)   |
| Application autotunes   | Simple autotune for BFB and DFTPP  |
| Data Analysis Software  | MassHunter and classic MSD ChemStation   |
| Application reports   | Environmental, drugs-of-abuse, aromatics in gasoline   |
| Custom reporting  | MassHunter: Microsoft Excel Classic MSD ChemStation: Custom Reporting  |
| Spectral libraries (optional)                                       | NIST, Wiley/NIST, Maurer/Pfleger/Weber   |
| Agilent Retention Time Locked<br>Mass Spectral Libraries (optional) | Pesticides and endocrine disrupter (including Japan Positive List), hazardous chemicals, indoor air toxics, forensic toxicology, metabolmics (Fiehn method), semi-environment volatiles, Solvents plus |
| Security  | Data and result security and tamper detection  |
| 21CFR11 Compliance  | MassHunter Quantitative Analysis and MassHunter GC/MS<br>Acquisition tools   |
| Multivariate analysis<br>(Optional)                                 | Mass Profiler Professional   |
| Sample class prediction (Optional)                                  | Sample Class Predictor or Mass Profiler Professional   |
| Other capabilities (optional)                                       | Deconvolution Reporting Software linked with RTL database for classic MSD ChemStation  |
|   |  |

### **Installation Checkout Specifications**

Agilent verifies GC/MSD system performance at the customer site. Test specifications are based on using Splitless Injection and an Agilent HP-5MSUI 30 m  $\times$  0.25 mm, 0.25 µm column, using helium as the carrier gas. Instrument Detection Limit (IDL) is a more effective metric of total system performance than signal-to-noise (S/N). Verification is an extensive (eight injections versus one) and reliable test to demonstrate system performance. See more about this type of test at http://www.chem.agilent.com/Library/technicaloverviews/Public/5990-8341EN.pdf

| 10 fg or less for Inert Extractor El source, Turbo Pump   |  |
|---|--|
| 24 fg or less for Inert El source, Turbo Pump   |  |
| <b>30 fg or less</b> for Inert El source, Diffusion Pump  |  |
| 8 sequential 1 $\mu L$ injections of 100 fg/ $\mu L$ 0FN $^4$ standard solution at nominal 272 u ion.                 |  |
| <b>1500:1 or higher</b> for Inert Extractor El source, Turbo Pump   |  |
| 600:1 or higher for Inert El source, Turbo Pump   |  |
| 300:1 or higher for Inert El source, Diffusion Pump   |  |
| 1- $\mu L$ injection of 1 pg/ $\mu L$ OFN standard scanning from 50 to 300 u at nominal 272 u ion.                    |  |
| 125:1 or higher   |  |
| 1-µL injection of 100-pg/µL BZP $^5$ standard scanning from 80 to 230 u at nominal 183 u ion                          |  |
| 600:1 or higher   |  |
| $2\text{-}\mu\text{L}$ injection of 100 fg/ $\mu\text{L}$ OFN standard scanning from 50 to 300 u at nominal 272 u ion |  |
|   |  |

<sup>&</sup>lt;sup>4</sup>Octafluoronaphthalene (OFN)

<sup>&</sup>lt;sup>5</sup>Benzophenone (BZP)

### Other Performance Specifications<sup>1</sup>

| PCI SIM S/N  | 10:1 or higher   |
|--|--|
| (Methane – reagent gas,  |  |
| Helium/Hydrogen carrier gas)   | $1\text{-}\mu\text{L}$ injection of 20-fg/ $\mu\text{L}$ BZP standard at nominal 183 u ion.                          |
| NCI SIM S/N  | 10:1 or higher   |
| (Methane, Helium/Hydrogen  |  |
| carrier gas)   | 1- $\mu L$ injection of 1-fg/ $\mu L$ OFN standard at nominal 272 u ion.   |
| PCI scan S/N (Ammonia)   | 500:1 or higher  |
|  | 1- $\mu L$ injection of 100-pg/ $\mu L$ BZP standard scanning from 80 to 230 u at nominal 183 u ion.                 |
| NCI scan S/N (Ammonia)   | 300:1 or higher  |
|  | $2\text{-}\mu\text{L}$ injection of 1-pg/ $\mu\text{L}$ OFN standard scanning from 50 to 300 u at nominal 272 u ion. |
| Both Extractor and Inert<br>EI scan S/N (H <sub>2</sub> carrier gas)<br>All EI sources | 100:1 or higher Turbo Pump<br>50:1 or higher Diffusion Pump  |
|  | $1\text{-}\mu L$ injection of $1\text{-}pg/\mu L$ OFN standard scanning from 50 to 300 u at nominal 272 u ion.       |

### **Ease-of-Maintenance**

- The GC inlet liner can be replaced in less than one minute without the use of tools with the Agilent 7890B GC system turn-top inlet.<sup>2</sup>
- A glass window shows ion source type, filament operation, and ion source connections.
- The source (including filaments and lenses) and electron multiplier can be removed from the instrument in less than one minute after venting.
- The optional micro ion gauge<sup>3</sup> can be replaced without removal of the mass spectrometer covers.

<sup>&</sup>lt;sup>1</sup>These specifications are reference of the performance. They are not confirmed at installation or familiarization.

<sup>&</sup>lt;sup>2</sup>Inlet temperature should be cool enough to touch when performing maintenance.

 $<sup>^3</sup>$ Micro ion gauge is shipped as standard for CI system but is available for EI systems as an option.

# Safety, Regulatory Compliance, and Operational Conditions

This instrument is manufactured and distributed under a quality system consistent with ISO13485. It complies with international regulatory, safety, and electromagnetic compatibility requirements. In addition, further testing was performed according to Agilent standards, to ensure operation after delivery and long-term usage.

| Safety                        | Canadian Standards Association (CSA): CAN/CSA-C22.2<br>No. 61010-1-04   |
|-------------------------------|---|
|                               | CSA/Nationally Recognized Test Laboratory (NRTL): UL 61010-1<br>International Electrotechnical Commission (IEC): 61010-1          |
|                               | EuroNorm (EN): 61010-1  |
| Electromagnetic compatibility | CISPR11/EN55011:Group1, Class A   |
| Sound emission                | EN 27779:1991 - sound pressure Lp <70 db  |
| Power                         | 120 V, 50/60 Hz, 1,100 VA; 200 V, 50/60 Hz, 1,100 VA;<br>220–240 V, 50/60 Hz, 1,100 VA  |
| Operating environment         | 15–35 °C, 40–80% relative humidity – noncondensing (operational) –40 °C–+70 °C, 0–95% relative humidity – noncondensing (storage) |

### Physical Requirements (with the Agilent 7890B GC)

| Dimensions (GC/MS) | 88 cm (w) $\times$ 56 cm(d) $\times$ 50 cm (h)<br>Additional space should be added for the auto injector, sample<br>tray, data system, and printer. Approximately 30 cm should be<br>available to the left of the MSD for ease of access to the system<br>in times of maintenance or repair. |
|--------------------|--|
| Weight (GC/MS)     | 100 to 112 kg not including automatic sampling systems (depending on configuration)  |

### **For More Information**

See the 5977A website at http://www/agilent.com/chem/5977A for further information.

This information subject to change without notice.

© Agilent Technologies, Inc., 2013 Published in USA, February 21, 2013 5991-1837EN

